



HPM-100-06/07

Ultra-High Speed Hybrid Detectors for TCSPC

Ultra fast instrument response function: <20 ps FWHM with SPC-180NX

HPM-100-06: 220 to 600 nm (Bialkali)

HPM-100-07: 220 to 850 nm (Multialkali)

No afterpulsing background

Excellent dynamic range of TCSPC measurements

Internal generators for PMT operating voltages

Power supply and control via bh DCC-100 card





Overload shutdown

Direct interfacing to all bh TCSPC systems



The HPM-100 module combines a Hamamatsu R10467 hybrid detector tube with a preamplifier and the generators for the tube operating voltages in one compact housing. The principle of the hybrid detector yields excellent timing resolution, a clean TCSPC instrument response function, high detection quantum efficiency, and extremely low afterpulsing probability. The absence of afterpulsing results in a substantially increased dynamic range of TCSPC measurements.

The HPM-100 module is operated via the bh DCC-100 detector controller of the bh TCSPC systems. The DCC-100 provides for power supply, gain control, and overload shutdown. The HPM-100 interfaces directly to all bh SPC or Simple Tau TCSPC systems. It is available with standard C-mount adapters, adapters for the bh DCS-120 confocal scanning FLIM system, and adapters for the NDD and BIG ports of the Zeiss LSM 710/780/880 NLO multiphoton laser scanning microscopes.

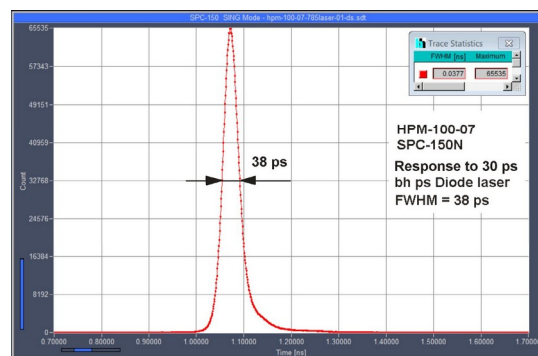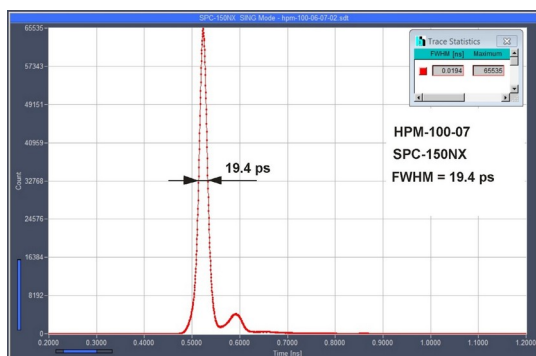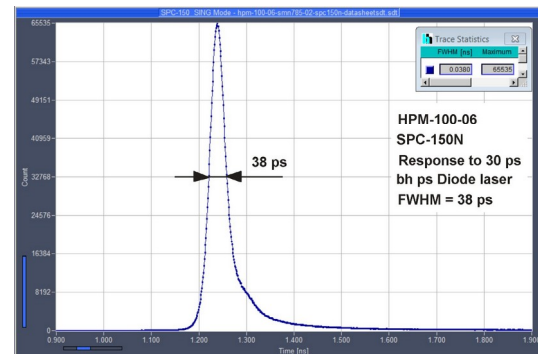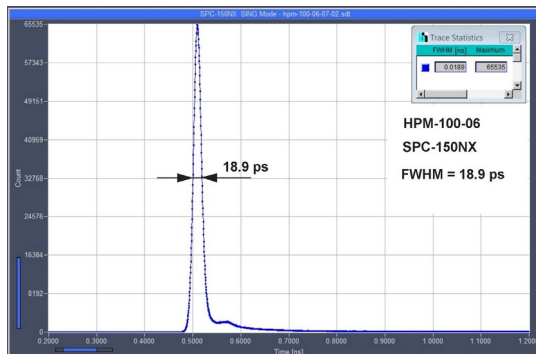


Left: Instrument response function, measured with 100-fs fibre laser. Recorded with SPC-150NX TCSPC module. Right: Response to pulses from bh picosecond diode laser, 30 ps pulse width. Recorded with SPC-150N TCSPC module.

Technology Leader in TCSPC



HPM-100-06/07

Detection quantum efficiency vs. wavelength

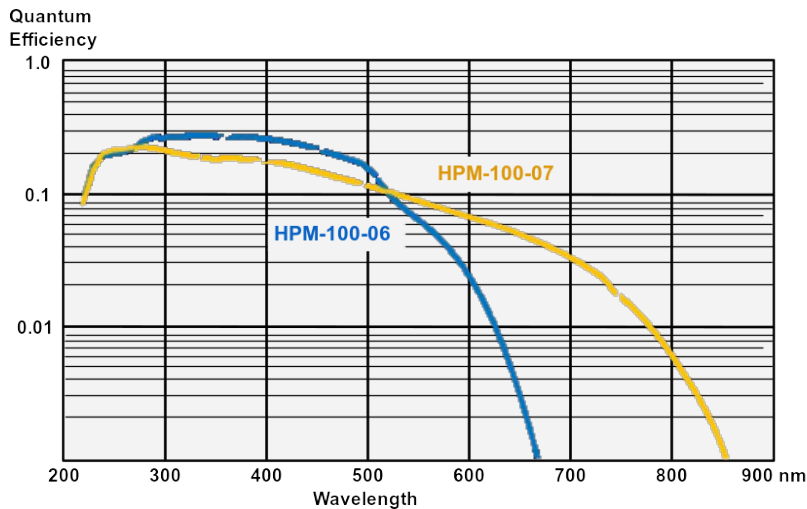

(after Hamamatsu Specifications)

Specifications, typical values

	-06 version	-07 version
Wavelength Range	220 nm to 650 nm	220 to 850 nm ¹⁾
Peak detection Quantum efficiency	28 % (at 350 nm)	16% at 400nm ¹⁾
Dark Count rate, T _{case} = 22°C	100 to 400 s ⁻¹	100 to 1000 s ⁻¹
Cathode Diameter	6 mm	3 mm
TCSPC IRF width (Transit Time Spread, with SPC-180NX)	<20 ps, FWHM	850 ps, FWHM
Single Electron Response Width	50 to 150 mV, -8000 V, V _{apd} 95% of V _{breakdown}	
Single Electron Response Amplitude		
Output Polarity		negative
Output Impedance		50 Ω
Max. Count Rate (Continuous)		10 MHz
Overload shutdown at		>15 MHz
Detector Signal Output Connector		SMA
Power Supply (from DCC-100 Card)		+ 12 V, +5 V, -12V
Dimensions (width x height x depth)		60 mm x 90 mm x 170 mm
Optical Adapters		C-Mount, DCS-120, LSM 710/780/880/980 NDD and BIG ports

1) according to Hamamatsu specifications

Related products: HPM-100-40/42 GaAsP and HPM-100-50 GaAs hybrid detector modules
Literature: The bh TCSPC Handbook, 9th edition, Becker & Hickl GmbH. Printed copies or electronic version on www.becker-hickl.com
 Sub-20ps IRF Width from Hybrid Detectors and MCP-PMTs. Application note, available from www.becker-hickl.com




Becker & Hickl GmbH
 Nunsdorfer Ring 7-9
 12277 Berlin, Berlin
 Tel. +49 / 30 / 787 56 32
 Fax. +49 / 30 / 787 57 34
 email: info@becker-hickl.com
www.becker-hickl.com



グローバルにネットワークを広げ、最先端の科学をお客様に提供

本社: 〒134-0088 東京都江戸川区西葛西6-18-14T.ビル Tel. 03-3686-4711
 営業所: 〒532-0003 大阪府大阪市淀川区宮原4-1-46 新大阪北ビル Tel. 06-6393-7411
 URL: <https://www.tokyoinst.co.jp> Mail: sales@tokyoinst.co.jp

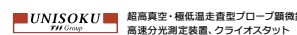

※本カタログに記載されている内容は、改良のため予告無く変更する場合があります。(製品の仕様、性能、価格などはカタログ発行当時のものです)
 ※本カタログに記載されている内容の一部または全部を無断で転載することは禁止されております。
 ※本カタログに記載されているメーカー名、製品名などは各社の商標または登録商標です。

International Sales Representatives



US:
Boston Electronics Corp
tcspc@boselec.com
www.boselec.com



UK:
Photonic Solutions PLC
sales@psplc.com
www.psplc.com



Japan:
Tokyo Instruments Inc.
sales@tokyoinst.co.jp
www.tokyoinst.co.jp



China:
DynaSense Photonics Co. Ltd.
info@dyna-sense.com
www.dyna-sense.com